Manual de uso

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www.medicallab.coreco.es

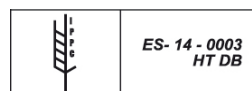
User's Manual

Refrigeration and freezing cabinets for **LABORATORIES** and **PHARMACIES**

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EN 63000:2018

EN ISO 22041:2019/A1:2019

Fabricante / *Manufacturer's name / Fabricant:*

CORECO S.A.

VAT ID: A14071559

Dirección / *Manufacturer's address / Adresse:*

CTRA. CÓRDOBA-MÁLAGA Km 80.800 - LUCENA (CÓRDOBA), SPAIN

Tipo de equipo / *Type of equipment / Type d'équipement:*

CABINETS FOR LABORATORIES AND PHARMACIES

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USER'S AND MAINTENANCE MANUAL

1. WARNING
2. GUARANTEE
3. INSTALLATION AND ASSEMBLY
4. TECHNICAL CHARACTERISTICS AND CONDITIONS OF USE
5. START-UP
6. RECOMMENDATIONS
7. PRECAUTIONS
8. MAINTENANCE
9. PERFORMANCE TESTS
10. AFTER-SALES FEEDBACK QUESTIONNAIRE
11. DECOMMISSIONING

INITIAL NOTICE

Before starting up the cabinet, we recommend reading and following the steps described in the instructions.

We would like to thank you for trusting us in purchasing this refrigerated unit. This manual is intended to ensure user safety and correct use and maintenance of the refrigeration unit. Please read it carefully before starting up your equipment. This refrigeration unit strictly complies with all applicable European directives and standards.

■ 1. WARNING

KEEP THIS MANUAL IN A SAFE PLACE WHERE YOU CAN REFER TO IT EASILY. IF IT BECOMES LOST, REQUEST A DUPLICATE, STATING THE MODEL, SERIAL NUMBER AND DATE OF PURCHASE.

THE MANUFACTURER ASSUMES THAT BOTH THE INSTALLATION TECHNICIAN AND THE USER OF THE UNIT HAVE THE BASIC TRAINING NECESSARY TO UNDERSTAND THE CONTENTS OF THIS MANUAL AND UNDERSTAND THE MAIN HYGIENE AND SAFETY STANDARDS.

REPAIR OPERATIONS MUST ONLY BE CARRIED OUT BY QUALIFIED TECHNICAL PERSONNEL.

ATTENTION: ONLY QUALIFIED PERSONNEL ARE AUTHORISED TO ACCESS THE ELECTRICAL PARTS OF THE CABINET, EITHER FOR INSTALLATION OR MAINTENANCE.

THESE PREREQUISITES SHALL BE AT THE USER'S EXPENSE:

- **PREPARATION OF THE PREMISES TO SUIT THE CABINET.**
- **MAINS POWER SUPPLY COMPLIANT WITH APPLICABLE REGULATIONS, ADEQUATE FOR THE CONSUMPTION AND SAFETY REQUIREMENTS OF THE UNIT.**
- **MATERIAL NECESSARY FOR THE CLEANING THE CABINET.**
- **WATER DRAINAGE POINT, IF NECESSARY FOR INSTALLATION OF THE REFRIGERATOR.**

THE MANUFACTURER DECLINES ALL RESPONSIBILITY IN THE FOLLOWING CASES:

- **IMPROPER USE OF THE CABINET.**
- **DAMAGE TO PERSONS OR OBJECTS DUE TO INCORRECT INSTALLATION.**
- **INSTALLATION NOT CARRIED OUT FOLLOWING THE PROCEDURES DESCRIBED IN THIS MANUAL.**
- **DEFECTS IN THE POWER SUPPLY.**
- **UNAUTHORISED MODIFICATIONS OR INTERVENTIONS.**
- **USE OF SPARE PARTS NOT SPECIFIC TO THE MODEL.**
- **TOTAL OR PARTIAL NON-COMPLIANCE WITH THE INSTRUCTIONS.**
- **INCIDENTS CAUSED BY THE TRANSPORT COMPANY OR WHEN MOVING THE UNIT AT THE INSTALLATION SITE.**

FOR ANY QUESTIONS, PLEASE CONTACT THE DISTRIBUTOR.

OTHER WARNINGS:

- **KEEP ALL VENTILATION OPENINGS IN THE APPLIANCE ENCLOSURE AND ANY HOUSING FREE OF OBSTRUCTIONS.**
- **DO NOT USE MECHANICAL DEVICES OR OTHER MEANS TO SPEED UP THE DEFROSTING PROCESS OTHER THAN THOSE RECOMMENDED BY THE MANUFACTURER.**
- **DO NOT DAMAGE THE COOLANT CIRCUIT.**
- **DO NOT USE ANY ELECTRICAL APPLIANCES INSIDE THE FOOD STORAGE COMPARTMENTS EXCEPT TYPES RECOMMENDED BY THE MANUFACTURER.**

■ 2. GUARANTEE

The device and its component parts are guaranteed one year from the date of sale. The guarantee covers replacement (freight prepaid) of the defective part or parts, provided that the defect is not due to misuse.

The guarantee does not cover labour for replacing parts.

The guarantee does not cover, even within the guarantee period, breakdowns or breakage of parts due to lack of maintenance or operation of the cabinet in working conditions other than those established in this manual, and the characteristics for each model listed in the catalogue (including power failures, high ambient temperatures, glass, lighting).

Returns **MUST BE AUTHORISED IN ADVANCE** by the manufacturer and must include the original packaging or similar.

The manufacturer's guarantee excludes any losses resulting from component failure.

■ 3. INSTALLATION AND ASSEMBLY

The cabinet should not be tipped up or laid down. If necessary, some appliances may be tipped up on the side indicated on the packaging but never on the other side.

It is advisable to unpack it on a firm, flat, stable floor. To move it to the installation site, use a wheelbarrow or pallet truck. Do not drag it across the floor. Ensure that the floor is not uneven.

Once it is in position, remove the packaging, leaving the unit free of all coverings and wrappings, avoiding knocks to the equipment. Now you can level the unit by turning the feet, up and down.

The unit is equipped with a cleaning drain which you should connect to a mains drain or a manually-removable tray.

IMPORTANT

If your unit is fitted with a backup battery for power failures, the switch to turn on battery power supply management to the controller is located at the **top rear** of the cabinet.

When you unpack the unit, the switch is set to "0" (battery management off).

When you connect the unit to the power supply (230V 50Hz), switch it to position "1" (battery management on).

Whenever you temporarily disconnect the machine for more than 12 hours, switch it back to the "0" position. When reconnecting to the mains (230V 50Hz), switch it back to position "1".

The following link takes you to information about each of the models: <https://medicallab.coreco.es/catalogo>

INITIAL CLEANING

First, remove the protective film from the steel (if present) using a non-sharp object. Clean off any adhesive residue with alcohol.

Absolutely avoid using tools or objects that could scratch or damage the unit. Once clean and dry, insert the accessories in the appropriate places.

We recommend cleaning the steel periodically with a sponge, hot water and neutral soap, then wiping it dry. **DO NOT USE DETERGENTS, PRODUCTS CONTAINING CHLORIDE OR SULPHUR OR ABRASIVE PRODUCTS.** Be especially careful not to get the electrical parts wet.

WARNING

When cleaning **STAINLESS STEEL** (if fitted), do not use iron or nickel scourers or even leave them resting on the surfaces, as they can leave ferrous residues that stick and cause rust due to contamination, compromising hygiene.

■ 4. TECHNICAL CHARACTERISTICS AND CONDITIONS OF USE

The cabinet's technical data, model and serial number are shown on the technical data plate provided on all models.

As well as its identifying information, the plate also states the unit's refrigerant gas and required electrical voltage.

		E.S. 14900-LUCENA(CORDOBA) Made in Spain (E.U.)		
N° SERIE SERIAL NR.				
MODELO Kode Mode	EQUIPO Equipement Equipment		AÑO CONSTRUCCIÓN Fabrikationsjahr Année de Fabrication Production Year	
CONDENSACION Kondensation Condensation		CONSUMO LÁMPARA Lampen Stromaufnahme Absorption lampe Lamp Absorption W		
POTENCIA FRIGORIFICA Gefriervermögen-Puissance-Frigorifique W Refrigeration Capacity		T Trabajo °C		
REFRIGERANTE Kältemittel-Refrigerant Type		g (± 10%)		
TENSION** Spannung-Tension V		Hz		
CONSUMO Stromaufnahme-Absorption Absorption		W max		
SISTEMAS CALEFACTORES W Heating Systems W Heizungsanlagen W Systèmes de Chauffage W		POTENCIA NOMINAL Nominleistung Puissance nominale W		
CLASE CLIMÁTICA Klima klasse Classe Climatique Climate Class		AGENTE EXPANSOR AISLAMIENTO Insulation expansion agent Isolation agent gonflant		

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The cabinet is designed for storing thermolabile products (substances subject to decomposition or change due to temperature) such as medicines, vaccines, reagents and physiological samples. When selecting the unit's temperature range, consider the following points:

- Type of product to be refrigerated or preserved.
- Room temperature.
- Frequency of door opening.
- Thermal insulation composition: 40 kg/m³ expanded injected polyurethane insulation with HFO blowing agent.

■ 5. START-UP

• INITIAL CLEANING

Clean the unit before connecting it to the mains. **(SEE INSTALLATION AND ASSEMBLY)**

• CHECKS

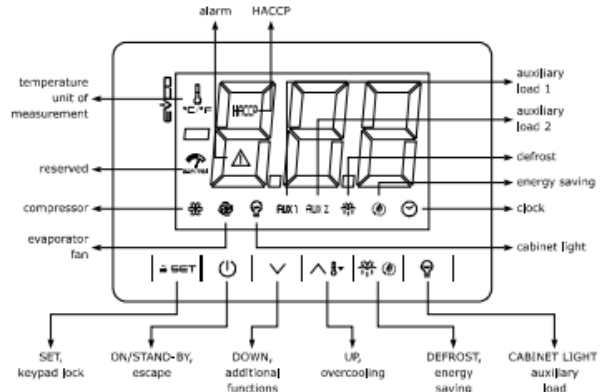
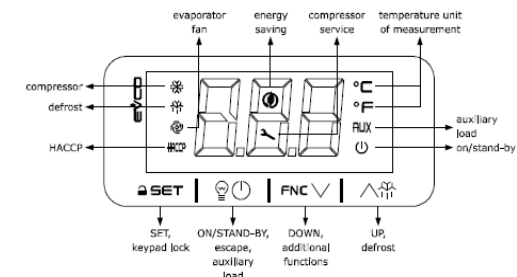
- Check that the mains voltage and frequency match those indicated on the unit's data plate.
- Never use the unit in corrosive or fire-hazard atmospheres.
- Check for any damage to the unit during transport. The manufacturer cannot be held liable for any damage during storage or caused by improper storage.
- Check the operation of the unit's moving parts.
- Check the state of the control panel, electrical wiring and connections.
- Check that the interior and adjoining fittings are present.

• GENERAL WIRING

- Check that the power cable is not damaged.
- If the power cable is damaged, replace it with another H05-VV-F or H05-VVH2-F type. This repair must be performed by qualified technical personnel.
- The mains power installation for the cabinet must have **EARTH GROUNDING**, and proper protection with a differential magneto-thermal switch.
- The cross-section of the mains power supply cable must be appropriate for the appliance's consumption.
- If several units are installed in line, they should be connected to the mains separately. Avoid using extension cables or multi-plug adaptors.
- The manufacturer declines all responsibility in case of non-observance of any of the above points.


- CONTROLLERS AND INDICATORS

DIGITAL CONTROLLERS

- Once the cabinet has been cleaned, connect it to the mains and press the on-off switch **(ON)**. The thermostat display indicates the air temperature inside the cabinet.
- Three minutes after switching it on the compressor LED should light up.
- The initial temperature shown on the thermostat display will be room temperature. As the machine continues operating, it will fall until it reaches the operating temperature.

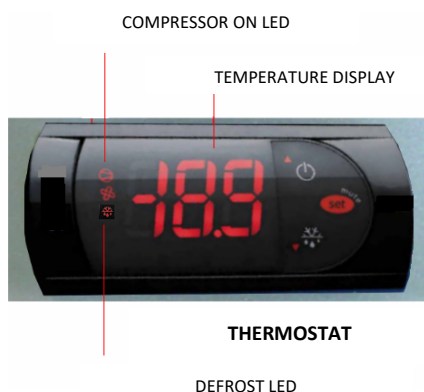SAFETY THERMOSTAT

On / Off

To turn on, press  until **ON** is displayed.

To turn off the unit, press the key again for 5 seconds until **OFF** is displayed, and the condenser unit will stop operating.

Set point temperature display.

Press  "set" for 1 second to display the set point. Release the button.

Press  the up arrow as necessary. Release the button.

Press  down arrow as necessary. Release the button.

To set the new value press . The ambient temperature probe reading is displayed.

Although you can select and reprogram the temperature using the thermostat, it is not advisable, as doing so may cause the unit to malfunction, altering the purpose for which it was programmed.

The unit user is liable for any damage caused by failure to comply with the observation described in the preceding point.

■ 6. RECOMMENDATIONS

- Position the product in such a way as to allow good air circulation, respecting the maximum load level shown inside the cabinet.
- Avoid opening the doors too often, as far as possible, and above all, never leave them open.

When the unit is to be turned off for a long time, the following steps are recommended:

- Remove products from inside the refrigerator.
- Turn off the main switch and unplug the unit.
- Clean the unit inside and outside.
- Leave the door ajar to allow air to circulate and avoid bad smells.

■ 7. PRECAUTIONS

- Although the unit is guaranteed to be stable, even with the doors open, never lean on the doors.
- Do not leave the cabinet outdoors.
- Do not alter the data plate or instructions provided by the manufacturer.
- Do not touch the equipment with wet or damp hands or feet.
- Do not touch the equipment while barefoot.
- Do not pull on the cable to unplug the unit.
- Avoid using adaptors, extension leads or multi-plug adaptors.
- Before performing any cleaning or maintenance operation, unplug the unit from the mains by first turning off the main switch and then the power plug.
- **NEVER** allow handling by children or persons with mental, sensory or intellectual disabilities. The same applies to persons who need supervision or instruction due to lack of experience or knowledge.
- Do not disassemble the protection of moving parts without first disconnecting them from the mains.
- Take the necessary precautions before putting your hands in the condenser compartment, as some parts may be hot, with the consequent risk of burns.

■ 8. MAINTENANCE

Before cleaning in any way, disconnect the cabinet from the mains power supply.

The unit has a drain for easier cleaning. When cleaning, you must remove the drain plug and clean the drain to prevent clogging with solids.

The proper functioning and durability of the unit depend largely on correct maintenance, carried out periodically.

The manufacturer disclaims any liability for non-compliance with the procedures described in this manual.

• GAS CONDENSER

Incorrect maintenance and not cleaning the gas condenser adequately can considerably reduce the refrigeration unit's performance and shorten the life of the compressor. The cleaner the condenser fins are, **the lower the energy consumption**.

This work should always be **done by a technical service**. If you are unsure, please contact your distributor. We recommend carrying out this operation at least every three months and at the beginning of each summer.

Before cleaning the gas condenser, unplug the unit and remove the protection grille or remove the condenser unit.

To remove dirt deposits on the condenser fins, any of the following procedures are recommended:

- Use a brush with soft bristles.
- Use a suitable vacuum cleaning device.
- Blow with compressed air to remove dirt.

- **EVAPORATORS**

Qualified technical personnel only

The only maintenance the evaporators require is adequate cleaning. This operation must only be done with the unit turned off. Wait for the ice between the fins to melt, or run a defrost cycle using the controller. This procedure can be done more quickly using warm water, but **NEVER** use hard or sharp objects.

To clean the battery or the casing, **DO NOT** use chemical degreasers with acid or bleach, chlorine or ammonia, or salts, as all these products aggressively attack copper and aluminium. Only use water with very dilute neutral detergents or products specially indicated for cleaning refrigeration equipment evaporators. The cleaner it is, **the greater the energy savings**.

- **AUTOMATIC EVAPORATION TRAY**

Qualified technical personnel only

Some units have an automatic evaporation tray for the water from defrosting the evaporator. The number and duration of refrigeration unit defrosts are factory set. The user must not change this setting.

- **GASKET**

The gasket provides a seal around the refrigerated cabinet doors and prevents the ingress of external elements.

A magnetic strip inside the rubber gasket holds it against the cabinet frame. Over time, grit particles may become stuck by the magnetic effect, which may oxidise, causing stains.

We recommend keeping the gasket free of foreign bodies that may produce rust, and of oily deposits, which over time may degrade the gasket rubber material.

■ 9. PERFORMANCE TESTS

- **CHECKS FOR MALFUNCTIONS**

Some malfunctions may have simple causes that the user can solve, so we recommend that you perform the following checks before calling in a specialised technician:

The device does not work:

- Check that it is plugged in correctly.
- Check that power is supplied to the socket.

The unit does not reach the cooling temperature:

- Check that there is no heat source nearby or that it is not exposed to direct sunlight.
- Check that the condenser grille is not obstructed.
- Check the set point temperature.
- Check that the product inside is correctly placed and does not block the air outlets of the internal fan. Check that the product has had enough time to cool.
- Check for ice on the evaporator.
- Check that the condenser is not clogged.
- Check that the unit is operating under normal conditions.
- Check that it is not exposed to drafts.

In case of strange or excessive noises:

- Check that the cabinet is levelled correctly; otherwise, it could vibrate.
- Check that no objects are rubbing against any of the refrigerator's moving parts.

- **NOISE TEST**

Carry out noise tests in a room with no noise-absorbing elements and away from large obstacles.

If the results of these tests are negative, please contact the service department in your area and provide the following information on the model, serial number and license plate number indicated on the data plate.

■ 10. AFTER-SALES FEEDBACK QUESTIONNAIRE

Scan the QR code or type the address in your browser's address bar:

<https://www.coreco.es/formularios/formulariopostventa>

■ 11. DECOMMISSIONING

Decommissioning

The refrigerated cabinet contains polyurethane foam, oil, plastic elements, metal parts and electrical and electronic components.

At the end of the cabinet's useful life, it must be collected for destruction or partial recovery. The user is responsible for delivering it to the collection centre specified by the local authority for professional WEEE recovery and recycling, always in compliance with applicable legal provisions.

The manufacturer is responsible for making the recovery, disposal and end-of-life treatment of the product feasible, either directly or through a collective system.

If the regulations are violated, specific sanctions may be applied, as established autonomously according to the legislation of each EC member state and binding on all those who are subject to such regulations.

Dismantling the Cabinet

While complying with each country's waste disposal regulations and fully respecting the natural environment, we recommend sorting the different cabinet parts by material so they can be disposed of separately or recovered.

This product's refrigerant circuit contains hydrocarbon refrigerant with a GWP value of 3 and does not contribute to global warming, so it can be released to the atmosphere.

Only release the refrigerant to the air outdoors.

RoHS Declaration of Conformity

Coreco hereby declares that the refrigerated cabinet referred to in this manual complies with the requirements of Directive 2002/95/EC/(RoHS).

In all homogeneous materials used in its manufacture, the possible presence of lead, mercury, hexavalent chromium, polybrominated biphenyls (PBB), and diphenyl ether (PBDE) is less than 0.01% by weight. This declaration is based on those by our suppliers of raw materials and components.

CORECO

Medical Lab

Ctr. Córdoba – Málaga, km 80'800

14.900 – Lucena (Córdoba) Spain

Tel.: +34 957 50 22 75

Fax: +34 957 51 42 98

info@coreco.es

www.medicallab.coreco.es